

Shim-pack Bio Diol & IEX Columns


Solution for Analysis of Peptides, Oligonucleotides and Biopharmaceuticals

The accurate analysis of biopharmaceuticals compounds is required for developing higher quality biopharmaceuticals. Shim-pack Bio Diol and IEX columns will improve the accuracy of the characterization of peptides, oligonucleotide and biopharmaceuticals.

Shim-pack Bio Diol : Size Exclusion Chromatography Columns

With different pore sizes, Shim-pack Bio Diol LC columns are effective for analysis of aggregates and fragments of mAb, oligonucleotides and carbohydrates.

Rapid mAb Aggregate Analysis using 2µm Shim-pack Diol-300 column

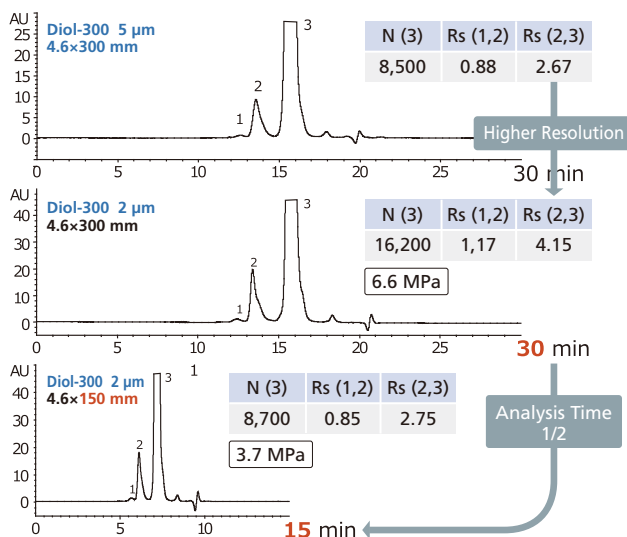

Shim-pack Bio Diol	Diol-60	Diol-120	Diol-200	Diol-300
Particle	Silica			
Ligand	Dihydroxypropyl(Diol)			
Particle Size	3 µm, 5 µm		2 µm, 3 µm, 5 µm	
Pore Size	6 nm	12 nm	20 nm	30 nm
pH Range	5.0 - 7.5			
Molecular Weight Range	below 10,000	1,000 - 100,000	5,000 - 300,000	20,000 - 1,000,000

By reducing the particle size from 5 µm to 2 µm, the resolution between aggregates and monomers was greatly improved. Furthermore, by reducing the column length from 300 mm to 150 mm using a 2 µm particle, 50% less run time was achieved, while maintaining resolution as compared to the original method with a 5 µm, 4.6x300 mm column.


Column	: Shim-pack Bio Diol-300
Eluent	: 0.1 M KH ₂ PO ₄ -K ₂ HPO ₄ (pH 7.0) with 0.2 M NaCl
Flow rate	: 0.2 mL/min
Column Temp.	: Ambient
Detection	: UV 280 nm
Sample	: Humanized monoclonal IgG1

Shim-pack Bio IEX : Ion Exchange Chromatography Columns

Shim-pack Bio IEX Columns are available in Q (quaternary ammonium) and SP (sulfopropyl) chemistries and are based on porous (Q and SP columns) and non-porous (Q-NP and SP-NP columns) hydrophilic polymers with low nonspecific adsorption. The porous particles offer excellent binding capacity with exceptionally high efficiency and the non-porous particles offer high efficiency and exceptional resolution.

Shim-pack Bio IEX	Q-NP	SP-NP	Q	SP
Particle	hydrophilic non-porous polymer		hydrophilic porous polymer	
Particle Size	3 µm, 5 µm		5 µm	
Ligand	- CH ₂ N ⁺ (CH ₃) ₃	- (CH ₂) ₃ SO ₃ ⁻	- CH ₂ N ⁺ (CH ₃) ₃	- (CH ₂) ₃ SO ₃ ⁻
pH Range	2 - 12			

Analysis of Synthesized Oligonucleotide (Single Strand DNA) using Shim-pack BIO IEX Q-NP

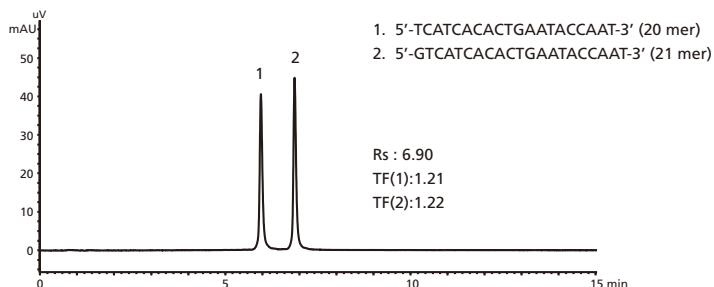

Column	: Shim-pack Bio IEX Q-NP 5 µm, 4.6x100 mm (P/N : 227-31003-03)
Mobile Phase A	: 10 mM NaOH
Mobile Phase B	: 10 mM NaOH with 1.0 M NaClO ₄
Gradient	: 25→55%B (0-15 min), 100%B (15-20 min)
Flow Rate	: 1.0 mL/min
Column Temp.	: 25 °C
Detection	: UV 260 nm
Inj. Volume	: 4 µL (5 nmol/mL)

Ordering Information

Shim-pack Bio Diol Columns

Particle Size	2 μ m		3 μ m			
Chemistry	Diol-200	Diol-300	Diol-60	Diol-120	Diol-200	Diol-300
4.6 \times 150 mm	227-31009-01	227-31010-01				
4.6 \times 300 mm	227-31009-02	227-31010-02	227-31007-01	227-31008-01	227-31009-03	227-31010-03
Max. Pressure	45 MPa		20 MPa			

Particle Size	5 μ m			
Chemistry	Diol-60	Diol-120	Diol-200	Diol-300
4.6 \times 300 mm	227-31007-02	227-31008-02	227-31009-04	227-31010-04
8.0 \times 300 mm	227-31007-03	227-31008-03	227-31009-05	227-31010-05
8.0 \times 30 mm (Guard Column)	227-31007-04	227-31008-04	227-31009-06	227-31010-06
Max. Pressure	20 MPa			
20 \times 300 mm	227-31097-01	227-31098-01	227-31099-01	227-31100-01
20 \times 500 mm	227-31097-02	227-31098-02	227-31099-02	227-31100-02
20 \times 50 mm (Guard Column)	227-31116-01	227-31117-01	227-31118-01	227-31119-01
Max. Pressure	10 MPa			

Shim-pack Bio IEX Columns

Shim-pack Bio IEX		Q-NP		SP-NP		Q	SP
Porality		Non-Porous				Porous	
Chemistry	Column Dimension	3 μm	5 μm	3 μm	5 μm	5 μm	
4.6×30 mm		227-31002-01	227-31003-01	227-31005-01	227-31006-01	227-31001-01	227-31004-01
4.6×50 mm		227-31002-02	227-31003-02	227-31005-02	227-31006-02	227-31001-02	227-31004-02
4.6×100 mm		227-31002-03	227-31003-03	227-31005-03	227-31006-03	227-31001-03	227-31004-03



For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

Company names, products/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation, its subsidiaries or its affiliates, whether or not they are used with trademark symbol "TM" or "®".

Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The contents of this publication are provided to you "as is" without warranty of any kind, and are subject to change without notice. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication.

Shimadzu Corporation

www.shimadzu.com/an/