

Application News

Analysis of Impurities in Alcohol-Based Hand Sanitizers by GC-MS

N. Iwasa

User Benefits

- ◆ This method can quantify impurities and alcohol (ethanol or isopropanol) in hand sanitizers.
- ◆ It is possible to measure listed impurities in a wide range of concentrations from low to high under a single analysis condition.
- ◆ FASST (Fast Automated Scan/SIM Type) analysis can identify compounds other than the listed impurities in hand sanitizers.

■ Introduction

Recently, various types of hand sanitizers have been produced to meet increasing demand due to the spread of infectious diseases, and the U.S. Food & Drug Administration (FDA) has announced a "Direct Injection Gas Chromatography Mass Spectrometry (GC-MS) Method for the Detection of Listed Impurities in Hand Sanitizers" (hereinafter, FDA hand sanitizer analysis method) for quality assessment of hand sanitizers. This method allows the evaluation of sanitizers with ethanol or isopropanol and can also be used in screening for "listed impurities" regulated under the FDA's "Guidance for Industry" on alcohol-based hand sanitizer products. Moreover, it is also possible to assay for % alcohol under the same analysis conditions as in the impurity analysis.

In this article, the quantitative analysis of impurities and alcohol concentration in ethanol-based sanitizers was conducted using a Shimadzu GCMS-QP2020 NX single quadrupole GC-MS, referring to the FDA hand sanitizer analysis method. The analysis conditions described in this article can be used to measure impurity and alcohol concentration in alcohol-based sanitizers. Furthermore, the listed impurities can be detected over a wide concentration range, and the system requirements were satisfied. It is also possible to identify compounds other than the listed impurities by a FASST analysis using high speed switching between the Scan mode and the SIM mode.

■ FDA Listed Impurities in Hand Sanitizers

The impurities listed in the FDA Guidance for hand sanitizer products are classified as Level 1 or Level 2, depending on the toxicity of the impurities in the sanitizer. The concentration of impurities should be below the FDA guideline limits. Table 1 shows the Level 1 and Level 2 impurities, their limit values and the concentration ranges for each of the measured target compounds which can be detected under the analysis conditions described in this article.

Table 1 FDA Listed Impurities and Detectable Concentration Ranges

| Compound Name | | Interim Limit Listed in FDA Guidances (ppm) | Concentration Ranges for this Method (µg/mL) |
|---------------|---------|---|--|
| Impurity | Level 1 | Methanol | NMT 630 |
| | | Benzene | NMT 2 |
| | | Acetaldehyde | NMT 50 |
| | | 1, 1-diethoxyethane | NMT 50 |
| | | Acetone | NMT 4400 |
| | Level 2 | 1-Propanol | NMT 1000 |
| | | Ethyl Acetate | NMT 2200 |
| | | 2-Butanol | NMT 6200 |
| | | Isobutanol | NMT 21700 |
| | | 1-Butanol | NMT 1000 |
| | | 3-Methyl-1-Butanol | NMT 4100 |
| | | Amyl Alcohol | NMT 4100 |
| | | Ethanol | - |
| | | Isopropyl Alcohol | - |

■ Analysis Conditions

The analytical conditions for GC and MS were listed in Table 2, 3 and 4.

Table 2 Instrument Configuration

| | |
|---------------|--|
| Model | : GCMS-QP2020 NX |
| Auto Injector | : AOC™-20i Plus |
| Auto Sampler | : AOC-20s Plus |
| Column | : SH-I-624Sil MS (30 m × 0.25 mm I.D., 1.4 µm) ^{*1} |

^{*1} P/N: 221-75962-30

Table 3 Analysis Conditions

| | |
|---------------------|--|
| GC | |
| Injection Temp. | : 250 °C |
| Injection Mode | : Split |
| Split Ratio | : 50 |
| Carrier Gas | : He |
| Carrier Gas Control | : Constant column flow rate (1 mL/min) |
| Column Temp. | : 40 °C (5 mins) - 30 °C/min - 240 °C (4 mins) ^{*1} |
| | Total 15.67 mins |
| Injection Volume | : 1.0 µL |
| MS | |
| Ion source Temp. | : 230 °C |
| Interface Temp. | : 240 °C |
| Ionization method | : EI |
| Measurement mode | : Scan/SIM (FASST mode) |
| Scan Range | : m/z 29 to 300 |
| SIM Ions | : Table 4 |
| Event time | : Scan 0.2 sec, SIM 0.3 sec |

^{*1} Under this oven heating condition, a high power oven model is necessary.

Table 4 MS Table

| Time (min) | Compound Name | Target | Ident 1 | Ident 2 |
|--------------|------------------------------|--------|---------|---------|
| 1.5 – 2.56 | Acetaldehyde | 44.0 | 29.0 | - |
| | Methanol | 31.0 | 29.0 | - |
| 2.56 – 4.00 | Ethanol | 31.0 | 45.0 | - |
| | Acetone | 43.0 | 58.0 | - |
| | Isopropyl Alcohol | 45.0 | 59.0 | - |
| 5.00 – 6.65 | 1-Propanol | 31.0 | 42.0 | 59.0 |
| | Ethyl Acetate | 43.0 | 61.0 | - |
| 6.65 – 7.20 | 2-Butanol | 45.0 | 59.0 | - |
| | Isobutanol | 43.0 | 42.0 | - |
| 7.20 – 10.00 | Benzene | 78.0 | 77.0 | - |
| | 1-Butanol | 56.0 | 41.0 | 31.0 |
| | 1, 1-diethoxyethane (Acetal) | 45.0 | 73.0 | 103.0 |
| | 3-Methyl-1-Butanol | 55.0 | 42.0 | 70.0 |
| | Amyl Alcohol | 42.0 | 55.0 | 41.0 |

The detector voltage was set to a relative value based on tuning.

■ Preparation of Standards and Samples

The FDA hand sanitizer analysis method describes methods for a spiked recovery test, impurities determination, and alcohol % assay.

The standards and samples were prepared in accordance with the FDA hand sanitizer analysis method. To represent the liquid and gel types of samples, the following three samples were employed in this study: two ethanol-based sanitizers, one commercially available off a store shelf (sanitizer ①) and the other produced by a sake brewer (sanitizer ②) both to represent the liquid form of sample and an ethanol-based sanitizer gel (sanitizer ③) to represent the gel form. The ethanol concentration of all of the sanitizer samples was approximately 80 % (v/v).

■ Spiked Recovery Test

The FDA method specifies a spiked recovery test for verification of the condition of the instruments and measurement samples. Therefore, measurement samples were prepared, and the analysis was carried out in specified injection order/number of measurements in accordance with the FDA hand sanitizer analysis method.

In evaluating the method precision, the system suitability criteria is that the %RSD of the peak area for each listed impurity for all injection (n=6) of the spiked recovery standard solution should be no more than 10 %. In evaluating the method's accuracy, the % Recovery for impurities should be within the range 80-120. The results of all measured target compounds satisfied these requirements.

As the results of the spiked recovery test, Table 5 shows the repeatability results of the spiked recovery standard solution and the spiked recovery test results for the three sanitizer samples.

Table 5 System Suitability and Spike and Recovery Test

| Compound Name | Peak area %RSD (n=6) | Recovery rate (%) | | |
|---------------------|-----------------------------------|-------------------|-------------|-------------|
| | Spiked recovery standard solution | Sanitizer ① | Sanitizer ② | Sanitizer ③ |
| Acetaldehyde | 1.633 | 96.3 | 91.8 | 107.9 |
| Methanol | 1.889 | 107.8 | 99.1 | 103.1 |
| Ethanol | 1.757 | - | - | - |
| Acetone | 1.737 | 107.2 | 102.9 | 103.8 |
| Isopropyl Alcohol | 1.755 | 109.7 | 104.8 | 105.7 |
| 1-Propanol | 1.802 | 105.7 | 107.1 | 100.3 |
| Ethyl Acetate | 2.035 | 107.3 | 101.9 | 100.8 |
| 2-Butanol | 1.956 | 104.3 | 102.6 | 101.7 |
| Isobutanol | 1.889 | 105.5 | 109.2 | 101.7 |
| Benzene | 8.014 | 106.0 | 90.2 | 104.7 |
| 1-Butanol | 2.255 | 105.7 | 103.3 | 100.4 |
| 1, 1-diethoxyethane | 1.700 | 110.9 | 104.4 | 103.7 |
| 3-Methyl-1-Butanol | 2.485 | 105.1 | 94.6 | 101.5 |
| Amyl Alcohol | 2.203 | 103.9 | 104.8 | 100.6 |

Due to the high concentration of ethanol in the sanitizer samples, the recovery rate was not calculated.

■ Impurities Determination

Determination of the impurities (Level 1 and Level 2) specified in the FDA Guidance was conducted. Measurement samples were prepared and the analysis was carried out by the specified injection order/number of measurements, in accordance with the FDA hand sanitizer analysis method.

As system suitability criteria, the %RSD of the peak area for each listed impurity for all injections of standard solution should be no more than 10 %. System suitability was satisfied for the all compounds subject to measurement.

The detailed analytical results of impurities determination are summarized in Table 6. The chromatogram of the standard solution and the SIM chromatograms of each compound are shown in Fig. 1 and 2 respectively.

Table 6 Results of Impurities Determination

| Compound Name | Standard solution | | Impurities in 100 mL of sanitizer (ppm) | | | FDA limit value NMT (ppm) |
|---------------------|-----------------------|----------------------|---|-------------|-------------|---------------------------|
| | Concentration (μg/mL) | Peak area %RSD (n=6) | Sanitizer ① | Sanitizer ② | Sanitizer ③ | |
| Acetaldehyde | 11.78 | 0.521 | 32.5 | 49.6 | ND | 50 |
| Methanol | 158.2 | 0.481 | ND | 31.4 | ND | 630 |
| Ethanol | 394.5 | 0.488 | - | - | - | - |
| Acetone | 158 | 0.290 | ND | ND | ND | 4400 |
| Isopropyl Alcohol | 392.5 | 0.265 | ND | ND | ND | - |
| 1-Propanol | 160.8 | 0.438 | ND | 827.8 | ND | 1000 |
| Ethyl Acetate | 180.4 | 0.537 | 16.0 | 168.1 | ND | 2200 |
| 2-Butanol | 161.6 | 0.549 | ND | ND | ND | 6200 |
| Isobutanol | 160.6 | 0.610 | ND | 648.6 | ND | 21700 |
| Benzene | 0.44 | 1.045 | ND | ND | ND | 2 |
| 1-Butanol | 162.0 | 0.432 | ND | 33.9 | ND | 1000 |
| 1, 1-diethoxyethane | 12.45 | 0.850 | 40.6 | 135.8 | ND | 50 |
| 3-Methyl-1-Butanol | 161.8 | 0.472 | ND | 1580.8 | ND | 4100 |
| Amyl Alcohol | 162.2 | 0.610 | ND | ND | ND | 4100 |

■ Alcohol Determination

Alcohol can also be assayed under the same analysis conditions as in the spiked recovery test and impurities determination. Measurement samples were prepared and the analysis was conducted by the specified injection order/number of measurements, in accordance with the FDA hand sanitizer analysis method.

Table 7 shows the %alcohol (v/v) results for ethanol in the three sanitizers.

* The FDA method states as follows: "If the ethanol or isopropanol peak for the impurity sample is more than 5x greater than the ethanol or isopropanol peak in the standard, dilute the impurity sample so that the resulting peak area should be approximately 0.5x the standard's peak area for those alcohols." In this article, the sanitizer samples were diluted 100 times with distilled water.

* Verification of system suitability criteria differs depending on whether verification is conducted on the same day as the impurity test or on a different day. For details, please refer to the FDA hand sanitizer analysis method.

Table 7 Determination of %Alcohol

| Quantitative value of ethanol (%) | | |
|-----------------------------------|-------------|-------------|
| Sanitizer ① | Sanitizer ② | Sanitizer ③ |
| 70.5 | 77.6 | 78.0 |

■ Calibration and Quantification

The FDA hand sanitizer analysis method includes the concentration range of each compound for this method shown in Table 1. Analysis conditions which enable analysis in this concentration range are necessary.

Six samples were prepared/measured by diluting the stock standard solution by 2, 5, 10, 20, 50, and 100 times. The linearity of calibration curves confirmed that measurement is possible over a wide range of concentrations.

Fig. 3 shows the calibration curves of each target compound.

* When saturating a high concentration sample or when the sensitivity of a low concentration sample is insufficient, adjust the detector voltage by approximately ± 0.1 .

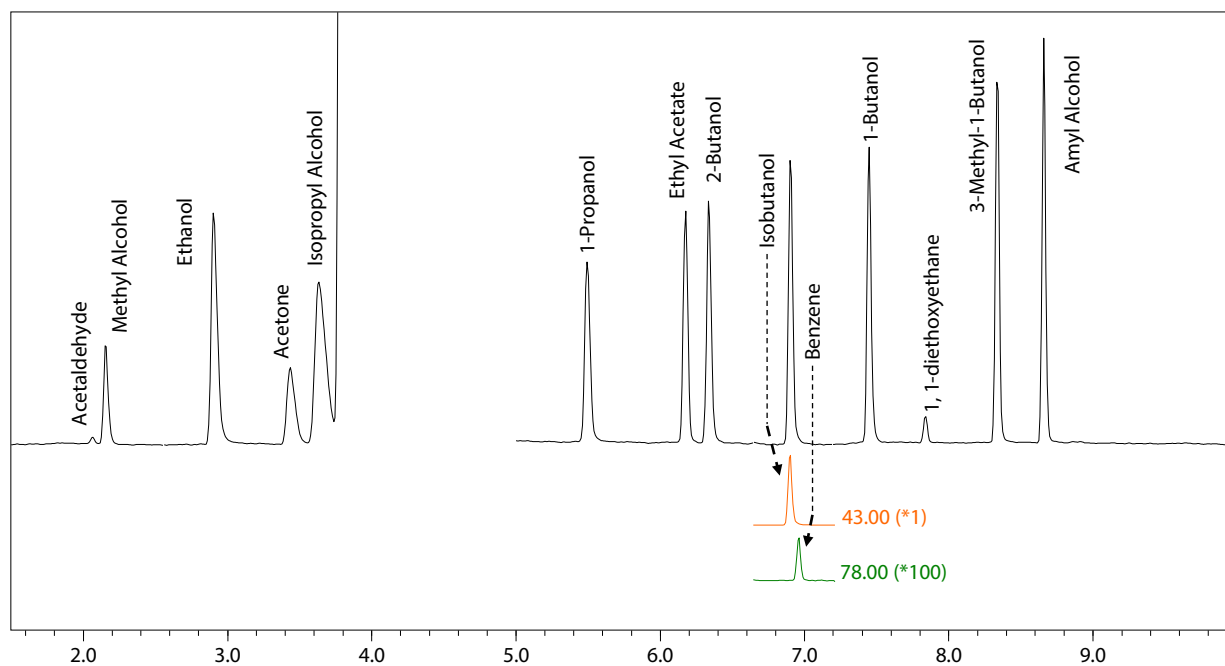


Fig. 1 TIC Chromatogram of Standard Solution (Isobutanol and Benzene: SIM Chromatogram)

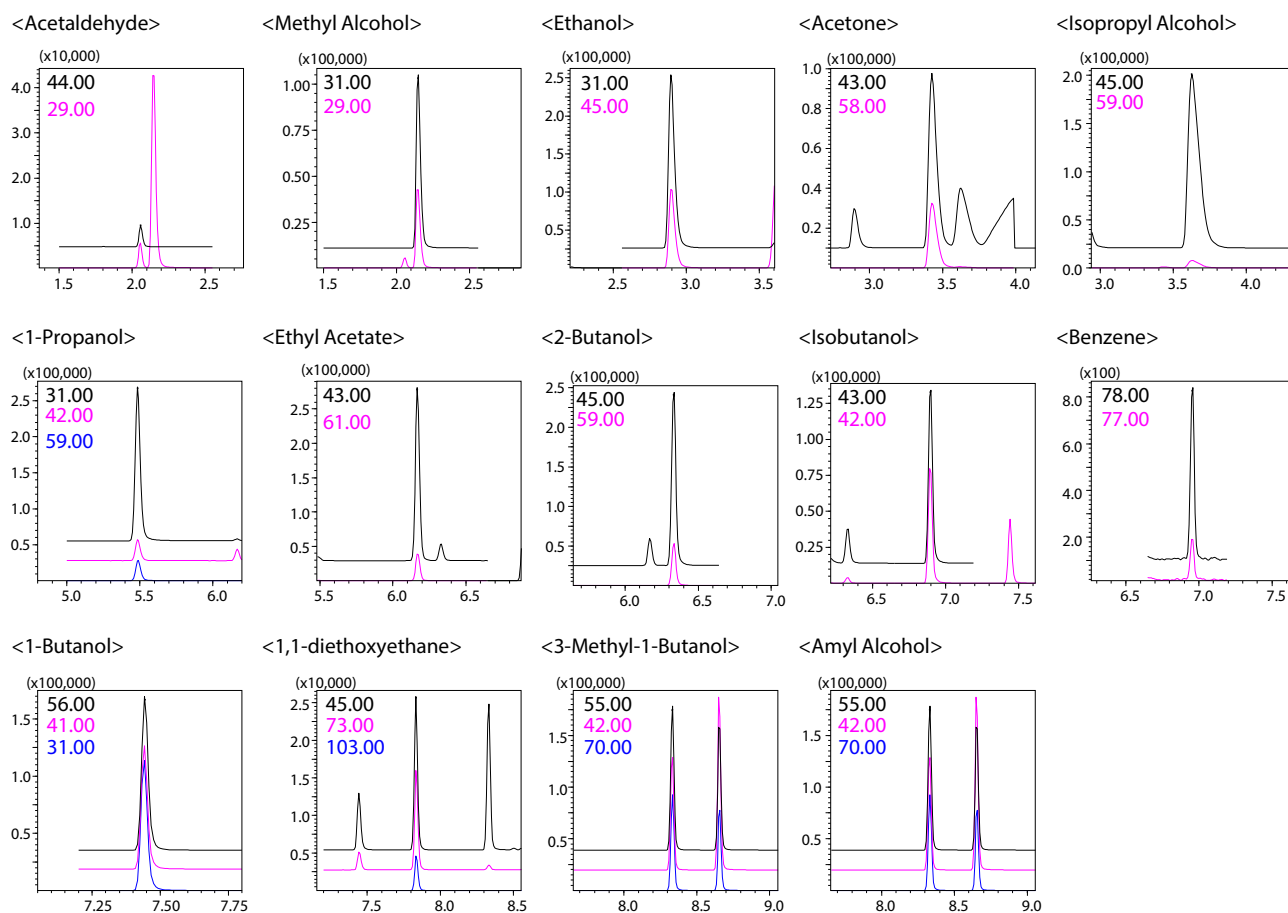


Fig. 2 SIM Chromatograms of Standard Solutions

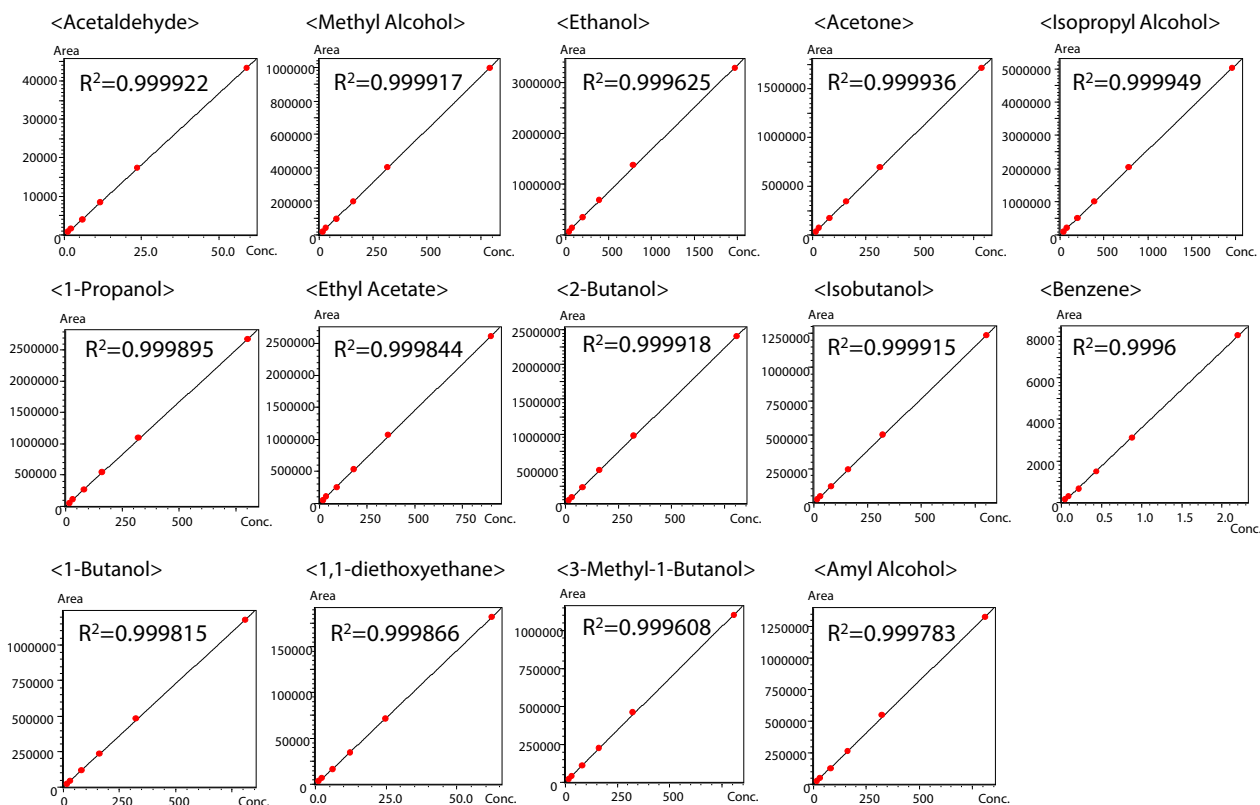


Fig. 3 Calibration Curve of Each Measured Compound

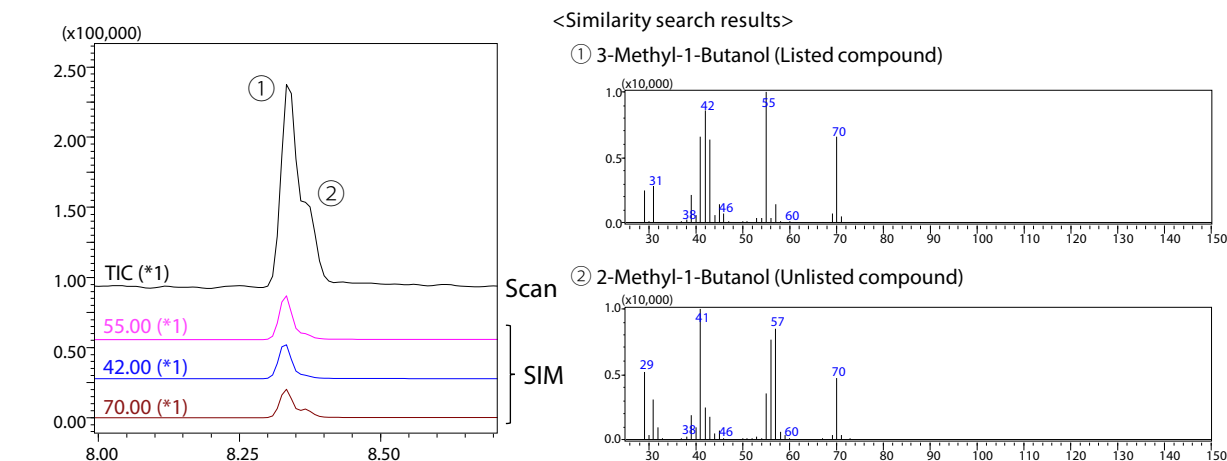


Fig. 4 Identification of Eluted Compounds other than Listed Impurities by FASST Mode

■ Identification by FASST

Using FASST (Fast Automated Scan/SIM Type) analysis, eluted compounds other than listed impurities can be identified. Among the sanitizer samples used in this experiment, a peak for a compound other than the listed impurities was detected in the sanitizer produced by a sake brewer (Sanitizer ②). This compound could be identified as 2-Methyl-1-Butanol from the Scan data. The results are shown in Fig. 4.

AOC is a trademark of Shimadzu Corporation or its affiliated companies in Japan and/or other countries.

■ Conclusion

GCMS-QP2020 NX can be used to quantify impurities and alcohol concentration in hand sanitizers. Identification of elution peaks not associated with listed impurities is also possible by FASST analysis.



Shimadzu Corporation

www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. See <http://www.shimadzu.com/about/trademarks/index.html> for details.

Third party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.